



Installation Instructions

TRILUTION® micro v2.4

Part Number 32000320

This document describes how to install TRILUTION® micro v2.4 on a PC.

RECOMMENDED PC CONFIGURATION

Processor	2.1 GHz 2-core
RAM	8.0 GB
Fixed Drive	400 GB
USB Port	One USB 2.0 or USB 3.0 port
Display	Resolution: 1440 x 900 (or greater) Color: True Color (32-bit) Text Size: 100%
Operating System	Microsoft® Windows® 7 Professional SP1 (64-bit) Microsoft® Windows® 10 Professional (64-bit)

Microsoft and Windows are registered trademarks of the Microsoft Corporation in the United States and/or other countries.

PRE-INSTALLATION CHECKLIST

Before beginning the installation, ensure that the PC name and domain are established. Changing these parameters after installation will result in errors.

- Ensure that the logged in user is a member of the Administrators Group.
- Close all running applications.
- Ensure that the format is set to match the same language and country as the operating system.
Note: After installation, to run TRILUTION micro, the Region Format needs to be set to English (United States).
- Temporarily disable antivirus software.
- Temporarily disable firewall.
- Temporarily disable checking for Windows Updates.
- Temporarily disable notifications.



INSTALLATION

One USB drive is used to install TRILUTION micro and the additional software that is required.

Note: If at any time during the installation there is a prompt to restart, then it is recommended to remove the USB drive before restarting.

The full installation will take approximately 30 minutes.

1. Connect the TRILUTION micro installation USB drive to a USB port on the PC.
2. Run **trilution_micro_pc_setup.exe** on the USB drive.
3. Follow the on-screen instructions to install any pre-requisites and then install TRILUTION micro.
 - While the installation is running, do not close any windows or boxes that appear. Take no action until prompted to.
 - If a User Account Control window appears during the installation, click **Yes**.
 - A restart may be required. After restarting, reconnect the USB drive. If the installation does not automatically resume, run **trilution_micro_pc_setup.exe**.
4. When the installation is finished, a prompt to restart the system appears. Disconnect the USB drive, and then restart the system. The system must be restarted before starting the software for the first time.

START TRILUTION® MICRO

Before starting TRILUTION micro for the first time, set the Region and Language Format to English (United States). Verify the following:

- The Decimal symbol used is a period and not a comma.
- The Short date and Long date formats are set to month/day/year.

The first time you start TRILUTION® micro, you must have local System Administrator privileges.

To start TRILUTION® micro:

- Windows® 7
 - Click the **Start** button and then select **All Programs > Gilson Applications > TRILUTION micro > TRILUTION micro x.x**.
- Windows® 10
 - Click the **Start** button and then select **(All apps) > Gilson Applications > TRILUTION micro x.x**.

If any Windows Security Alerts appear with Gilson, Inc. as the Publisher, click **Unblock**.

LOG IN

To log in to TRILUTION micro for the first time, enter the following User name and Password and then select **Accept**.

User name: admin

Password: Gilson268

The Password is case sensitive and must be entered exactly.

INSTALLATION QUALIFICATION

Complete the supplied Installation Qualification document.